

Алматы (7273)495-231
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Россия +7(495)268-04-70

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Казахстан +7(7172)727-132

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Нижний Новгород (831)429-08-12
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Киргизия +996(312)96-26-47

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Ульяновск (8422)24-23-59
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

<https://biochromato.nt-rt.ru> || boa@nt-rt.ru

- DMSO, DMF, WATER can be evaporated with NO BUMPING
- Solvent removal directly from HPLC / LC / GC vials
- Sample concentration after fractionation, NMR measurement, etc.

Smart Evaporator

MADE IN JAPAN

Over 1600 Units Sold Worldwide



In Vial Concentration




Grand Prize

Grand Prize

Award of Excellence

DMSO, DMF, Water

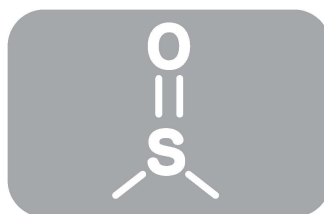

The 42nd Invention Award
(the Japan Society for the Advancement of Inventions)

Bump Free Technology



The 7th Kanagawa "Industrial Navi Award"
(Kanagawa Welfare Foundation for Entrepreneurs, Japan)

The 26th Small Business Excellence,
New Technology / New Product Award
(Resona Foundation / NK Industrial R.I., Japan)

In Vial Concentration




The Spiral Plugs come in various sizes to fit to container mouth inner diameter from 4mm to 32mm. With the aluminum beads on the bottom*, various sizes and shapes of tubes / vials are applicable for concentration.

Containers from other instruments or storage can be used directly, minimizing the loss by sample transfer (*features for C1 and K4)



DMSO, DMF, Water

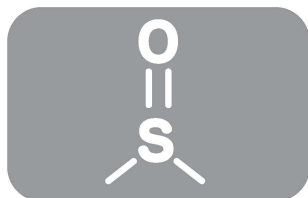

The Smart Evaporator enables efficient evaporation of high-boiling point solvents - even DMSO, DMF, and water.

All models are with the heating function, and the covered models* allow the use of inert gas, which accelerates faster concentration. (*C1 and C10)

Fast and Effective Evaporation

Solvent / Capacity	Flow rate	40°C	70°C
Hexane / 5 mL (20 mL vial)	15L/min	3 min	3 min
Acetone / 5 mL (20 mL vial)	15L/min	4 min	3 min
Ethyl acetate / 5 mL (20 mL vial)	15L/min	6 min	5 min
Acetonitrile / 5 mL (20 mL vial)	15L/min	9 min	8 min
Methanol / 5 mL (20 mL vial)	15L/min	10 min	6 min
Ethanol / 5 mL (20 mL vial)	15L/min	10 min	6 min
*Deionized Water / 5 mL (20 mL vial)	29~30L/min	49 min	28 min
*DMF / 5 mL (20 mL vial)	20~30L/min	43 min	24 min
*DMSO / 5 mL (20 mL vial)	23~30L/min	195 min	69 min
*NMP / 5 mL (20 mL vial)	25~30L/min	N/A	71 min

* Gas purge data


Bump Free Technology




The novel Spiral Plug technology generates a helical flow of air over the surface of the solvent, thereby increasing the surface area.

Since the container isn't under high vacuum, there is no risk of bumping or splashing.

Vacuum Vortex Concentration

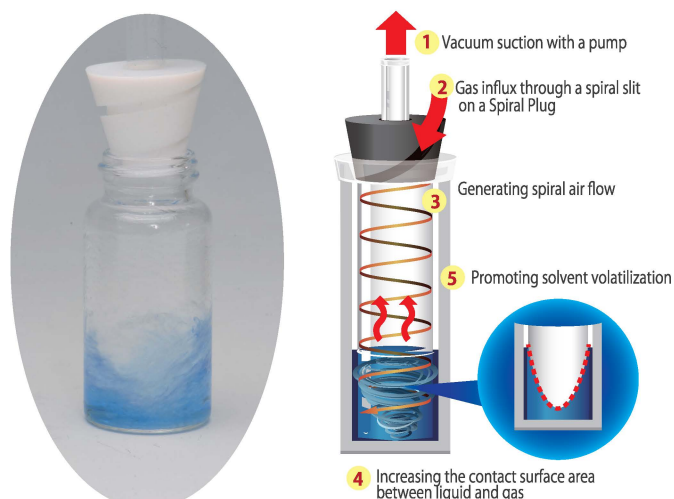


Successful Applications with the Smart Evaporator

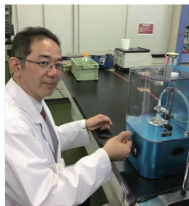

DMSO and DMF can be evaporated so easily and efficiently under 40°C - amazing!
I'm also happy with the efficiency in the sample recovery.

- Dr. Matsufuji, Food Analysis, Nihon University

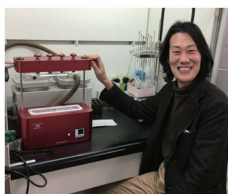

It is great being able to flexibly concentrate fractionated samples. No preparation time, each sample can be set when ready - we are saving a lot of time.

- Dr. Koyama, Food Chemistry and functional Nutrition, Tokyo University of Marine Science and Technology



Even the fraction container can be used for concentration.

The sample can be evaporated in the containers that we use, then directly injected into mice with a syringe.

We now plan the concentration during lunch break, and radioisotope with 2hours' half life can be efficiently tested.

- Dr. Hanaoka, Bioimaging Information Analysis, Gunma University



Users' Fields

Bioimaging, Biomaterial, Chemical Biology,
Chemical Bioscience, Molecular Biology,
Nuclear Medicine, Forensic Medicine, Drug Assay,
Organic Synthesis, Pharmacognosy,
Archaeometry, Natural Products Chemistry,
Nutritional Biochemistry, Food Analysis,
Catalyst Chemistry, Surface Chemistry,
Complex Chemistry, Inorganic, Chemistry,
Particle Chemistry, Material Chemistry,
Metalorganic Chemistry, Analytical Chemistry,

Application Examples

- Solvent removal (for substitution) after HPLC, LC/MS, GC/MS
- Sample concentration after preparative fraction
- Solvent removal after NMR measurement
- Concentration of natural product extract
- Concentration after solid phase extraction
- Concentration of radioisotope labeled compounds
- Concentration of surfactant solutions
- Pretreatment for multiple sample in drug analysis
- Pretreatment for analysis of photosynthetic pigment

Installation example

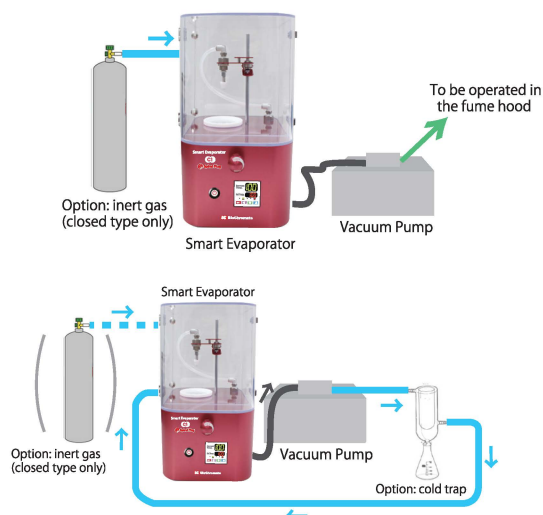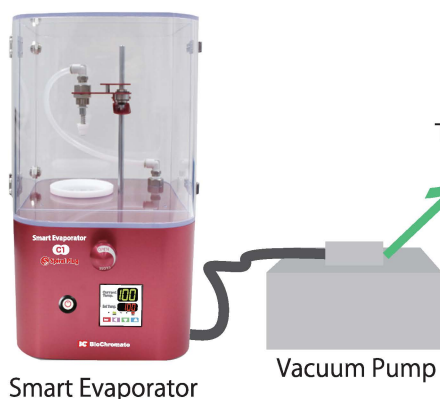

Smart Evaporator – models, accessories & specs

Smart Evaporator C1

Catalog code	CEV1B-SU / SQ / GR / SK
Size	W220 x D227 x H382 mm
Weight	4.3kg
Heating	Room temp. ~ 100C
Heating cup	Round shape / ϕ 69mm
Color	4 (Sun, Squall, Grass, Sky)
Gas purge	Yes
CE marking	Yes



Smart Evaporator K4

Smart Evaporator C10

Catalog code	CEK4-SU-P (plug size 1-5)	CEV10L-(color)-P (plug size 1-3)
Size	W434 x D156 x H410mm	W350 (480) x D300 x H531 mm
Weight	4.8kg	20kg
Heating	Room temp. ~ 100C	
Heating cup	Rectangle / W226 x D55mm	Round shape / with vial holder
Color	1 (Sun)	4 (Sun, Squall, Grass, Sky)
Gas purge	No (special order possible)	Yes
CE marking	No	Yes



Spiral Plug SP-S series

Spiral Plug SP-V series

Catalog code	SP-P (plug size No. 1-5)	SP-V (plug size No. 1-5)
Plug Material	PTFE	fluoroelastomer (FKM)
Size No. (Inner ϕ of container)	1 (4~7mm) 2 (7~11mm) 3 (11~17mm) 4 (15~24mm) 5 (24~32mm)	



Filter for C1

Catalog code	CEV-F-00
Size	L110 x ϕ 26 mm
Connection	R1/4 (male)
Filter material	Polyethylene



4Kit for C1

Catalog code	UW4-C1B
Parts Included	4-channel head 3-Way Vial Holder 4 x SP-P1 4 x SP-P2 4 x 2ml Vial 4 x 4ml Vial 4 x 6ml Vial



One kit converts C1 to 4 position that accommodates vials of 3 sizes

Pump Information

All Smart Evaporators need to be connected to a DIAPHRAGM VACUUM pump for operation*. The pump has to be capable of continuous suction at atmospheric pressure; and should reach the pressure of -40kPa (61.3×10^3 Pa) ~ -50kPa (51.3×10^3 Pa) or more.

The reference flow rate per plug size (= container mouth inner diameter) is as shown below. For multi-channel models, multiply the flow rate by the number of the channel.

*Please contact us for more detail and recommendations.

Inner diameter of container	Plug Size	Reference suction flow rate
4~7mm	Plug Size:1	13L/min
7~11mm	Plug Size:2	15L/min
11~17mm	Plug Size:3	30~33L/min
15~24mm	Plug Size:4	50L/min
24~32mm	Plug Size:5	58L/min

Алматы (7273)495-231	Иваново (4932)77-34-06	Магнитогорск (3519)55-03-13	Ростов-на-Дону (863)308-18-15	Тольятти (8482)63-91-07
Ангарск (3955)60-70-56	Ижевск (3412)26-03-58	Москва (495)268-04-70	Рязань (4912)46-61-64	Томск (3822)98-41-53
Архангельск (8182)63-90-72	Иркутск (395)279-98-46	Мурманск (8152)59-64-93	Самара (846)206-03-16	Тула (4872)33-79-87
Астрахань (8512)99-46-04	Казань (843)206-01-48	Набережные Челны (8552)20-53-41	Санкт-Петербург (812)309-46-40	Тюмень (3452)66-21-18
Барнаул (3852)73-04-60	Калининград (4012)72-03-81	Нижний Новгород (831)429-08-12	Саратов (845)249-38-78	Ульяновск (8422)24-23-59
Белгород (4722)40-23-64	Калуга (4842)92-23-67	Новокузнецк (3843)20-46-81	Севастополь (8692)22-31-93	Улан-Удэ (3012)59-97-51
Благовещенск (4162)22-76-07	Кемерово (3842)65-04-62	Ноябрьск (3496)41-32-12	Саранск (8342)22-96-24	Уфа (347)229-48-12
Брянск (4832)59-03-52	Киров (8332)68-02-04	Новосибирск (383)227-86-73	Симферополь (3652)67-13-56	Хабаровск (4212)92-98-04
Владивосток (423)249-28-31	Коломна (4966)23-41-49	Омск (3812)21-46-40	Смоленск (4812)29-41-54	Чебоксары (8352)28-53-07
Владикавказ (8672)28-90-48	Кострома (4942)77-07-48	Орел (4862)44-53-42	Сочи (862)225-72-31	Челябинск (351)202-03-61
Владимир (4922)49-43-18	Краснодар (861)203-40-90	Оренбург (3532)37-68-04	Ставрополь (8652)20-65-13	Череповец (8202)49-02-64
Волгоград (844)278-03-48	Красноярск (391)204-63-61	Пенза (8412)22-31-16	Сургут (3462)77-98-35	Чита (3022)38-34-83
Вологда (8172)26-41-59	Курск (4712)77-13-04	Петрозаводск (8142)55-98-37	Сыктывкар (8212)25-95-17	Якутск (4112)23-90-97
Воронеж (473)204-51-73	Курган (3522)50-90-47	Псков (8112)59-10-37	Тамбов (4752)50-40-97	Ярославль (4852)69-52-93
Екатеринбург (343)384-55-89	Липецк (4742)52-20-81	Пермь (342)205-81-47	Тверь (4822)63-31-35	
Россия +7(495)268-04-70		Казахстан +7(7172)727-132	Киргизия +996(312)96-26-47	

https://biochromato.nt-rt.ru || boa@nt-rt.ru